

Stock Solution



TD-S Revision 2.0

Creation Date: 4/8/2019

Revision Date: 4/8/2019

Coelenterazine Stock Solution for *In Vitro* Use

Instructions

1. Add 1 mg coelenterazine ([GoldBio Catalog # CZ](#)) to 1 ml of acidified methanol (add 50 μ l of concentrated HCl to 10 ml methanol).
2. Aliquot and store at -80°C for up to 4 weeks.
3. For use, thaw tube and add 25.41 μ l of coelenterazine stock solution to 40 ml DPBS (1.5 μ M final concentration).

Gold Biotechnology

St. Louis, MO

Ph: (800) 248-7609

Web: www.goldbio.com

Email: contactgoldbio86@goldbio.com